

Flow Cell Wash Kit (EXP-WSH003)

Version: WFC_9088_v1_revE_18Sep2019
 Last update: 12/03/2020



Flow Cell Number:

DNA Samples:

Before start checklist		
Materials	Consumables	Equipment
<input type="checkbox"/> Flow Cell Wash Kit (EXP-WSH003)		<input type="checkbox"/> Ice bucket with ice
		<input type="checkbox"/> Pipettes and pipette tips P20, P1000
INSTRUCTIONS		NOTES/OBSERVATIONS
Flushing a MinION/GridION Flow Cell		
<p>Preparation to run the washing procedure.</p> <ul style="list-style-type: none"> <input type="checkbox"/> Place the tube of Wash Solution A on ice. Do not vortex the tube. <input type="checkbox"/> Thaw one tube of Wash Solution B at RT. <input type="checkbox"/> Mix the contents of Wash Solution B thoroughly by vortexing, spin down briefly and place on ice. <p>In a clean 1.5 ml Eppendorf DNA LoBind tube, prepare the following Wash Mix:</p> <ul style="list-style-type: none"> <input type="checkbox"/> 20 µl Wash Solution A (A) <input type="checkbox"/> 380 µl Wash Solution B (B) <ul style="list-style-type: none"> <input type="checkbox"/> Mix well by pipetting, and place on ice. Do not vortex the tube. <input type="checkbox"/> Stop or pause the sequencing experiment in MinKNOW, and leave the flow cell in the device. <input type="checkbox"/> Ensure that the priming port cover and SpotON sample port cover are in the positions indicated in the figure below. <input type="checkbox"/> Using a P1000, remove all fluid from the waste channel through Waste port 1. As both the priming port and SpotON sample port are closed, no fluid should leave the sensor array area. 		
<p>IMPORTANT</p> <ul style="list-style-type: none"> <input type="checkbox"/> It is vital that the flow cell priming port and SpotON sample port are closed to prevent air from being drawn across the sensor array area, which would lead to a significant loss of sequencing channels. 		
<ul style="list-style-type: none"> <input type="checkbox"/> Rotate the flow cell priming port cover clockwise so that the priming port is visible. <p>Check for air between the priming port and the sensor array. If necessary, using a P1000 draw back a small volume to remove any air (a few µls):</p> <ul style="list-style-type: none"> <input type="checkbox"/> Set a P1000 pipette to 200 µl <input type="checkbox"/> Insert the tip into the priming port <input type="checkbox"/> Turn the wheel until the dial shows 220-230 µl, or until you can see a small volume of buffer/liquid entering the pipette tip. <input type="checkbox"/> Visually check that there is continuous buffer from the priming port across the sensor array. 		
<p>IMPORTANT</p> <ul style="list-style-type: none"> <input type="checkbox"/> Take care when drawing back buffer from the flow cell. Do not remove more than 20-30 µls, and make sure that the array of pores are covered by buffer at all times. Introducing air bubbles into the array can irreversibly damage pores. 		

Flow Cell Wash Kit (EXP-WSH003)

Version: WFC_9088_v1_revE_18Sep2019
 Last update: 12/03/2020



Flow Cell Number:

DNA Samples:

INSTRUCTIONS	NOTES/OBSERVATIONS
<ul style="list-style-type: none"> <input type="checkbox"/> Load 400 µl of the prepared Wash Mix into the flow cell via the priming port, avoiding the introduction of air. <input type="checkbox"/> Close the priming port and wait for 30 minutes. <input type="checkbox"/> Ensure that the priming port cover and SpotON sample port cover are in the positions indicated in the figure below. <input type="checkbox"/> Using a P1000, remove all fluid from the waste channel through Waste port 1. As both the priming port and SpotON sample port are closed, no fluid should leave the sensor array area. 	
<p>IMPORTANT</p> <ul style="list-style-type: none"> <input type="checkbox"/> It is vital that the flow cell priming port and SpotON sample port are closed to prevent air from being drawn across the sensor array area, which would lead to a significant loss of sequencing channels. 	
<p>Follow one of the two options described in the next steps of the protocol.</p>	
<p>To run a second library on a MinION/GridION flow cell straight away</p>	
<p>IMPORTANT</p> <ul style="list-style-type: none"> <input type="checkbox"/> The buffers used in this process are incompatible with conducting a Flow Cell Check step prior to loading the subsequent library. In order to check your flow cell, follow the instructions in the next section "To store the MinION/GridION/PromethION flow cell for later use" before priming and loading the flow cell. <input type="checkbox"/> To run a second library straight away, follow the instructions in the "Priming and loading the flow cell" section of the relevant protocol. 	
<p>To store the MinION/GridION flow cell for later use</p>	
<ul style="list-style-type: none"> <input type="checkbox"/> Thaw one tube of Storage Buffer (S) at RT. <input type="checkbox"/> Mix contents thoroughly by pipetting and spin down briefly. <input type="checkbox"/> Rotate the flow cell priming port cover clockwise so that the priming port is visible. <p>Check for air between the priming port and the sensor array. If necessary, using a P1000 draw back a small volume to remove any air (a few µls):</p> <ul style="list-style-type: none"> <input type="checkbox"/> Set a P1000 pipette to 200 µl <input type="checkbox"/> Insert the tip into the priming port <input type="checkbox"/> Turn the wheel until the dial shows 220-230 µl, or until you can see a small volume of buffer/liquid entering the pipette tip. <input type="checkbox"/> Visually check that there is continuous buffer from the priming port across the sensor array. <ul style="list-style-type: none"> <input type="checkbox"/> Slowly add 500 µl of Storage Buffer (S) through the priming port of the flow cell. <input type="checkbox"/> Close the priming port. <input type="checkbox"/> Using a P1000, remove all fluid from the waste channel through Waste port 1. As both the priming port and SpotON sample port are closed, no fluid should leave the sensor array area. 	

Flow Cell Wash Kit (EXP-WSH003)

Version: WFC_9088_v1_revE_18Sep2019
 Last update: 12/03/2020



Flow Cell Number:

DNA Samples:

INSTRUCTIONS	NOTES/OBSERVATIONS
<p>IMPORTANT</p> <p><input type="checkbox"/> It is vital that the flow cell priming port and SpotON sample port are closed to prevent air from being drawn across the sensor array area, which would lead to a significant loss of sequencing channels.</p>	
<p><input type="checkbox"/> The flow cell can now be stored at 4-8° C.</p> <p><input type="checkbox"/> When you wish to reuse the flow cell, remove the flow cell from storage, and allow it to warm to RT for ~5 minutes.</p>	
<p>Flushing a PromethION flow cell</p>	
<p>Preparation to run the washing procedure.</p> <p><input type="checkbox"/> Place the tube of Wash Solution A on ice. Do not vortex the tube.</p> <p><input type="checkbox"/> Thaw one tube of Wash Solution B at RT.</p> <p><input type="checkbox"/> Mix the contents of Wash Solution B thoroughly by vortexing, spin down briefly and place on ice.</p> <p>In a clean 1.5 ml Eppendorf DNA LoBind tube, prepare the following Wash Mix:</p> <p><input type="checkbox"/> 20 µl Wash Solution A (A)</p> <p><input type="checkbox"/> 380 µl Wash Solution B (B)</p> <p><input type="checkbox"/> Mix well by pipetting, and place on ice. Do not vortex the tube.</p> <p><input type="checkbox"/> Stop or pause the sequencing experiment in MinKNOW, and leave the flow cell in the device.</p> <p><input type="checkbox"/> Rotate the inlet port cover clockwise to reveal the inlet port.</p> <p>A small tract of air may be visible beyond the inlet port. If necessary, using a P1000 draw back a small volume to remove any air (a few µls):</p> <p><input type="checkbox"/> Set a P1000 pipette to 200 µl</p> <p><input type="checkbox"/> Insert the tip into the inlet port</p> <p><input type="checkbox"/> Turn the wheel until the dial shows 220-230 µl, or until you can see a small volume of buffer/liquid entering the pipette tip.</p> <p><input type="checkbox"/> Load 400 µl of the prepared Wash Mix into the flow cell via the inlet port, avoiding the introduction of air.</p> <p><input type="checkbox"/> Close the inlet port and wait for 30 minutes.</p>	
<p>Follow one of the two options described in the next steps of the protocol.</p>	
<p>To run a second library on a PromethION flow cell straight away</p>	
<p>IMPORTANT</p> <p><input type="checkbox"/> The buffers used in this process are incompatible with conducting a Flow Cell Check step prior to loading the subsequent library. In order to check your flow cell, follow the instructions in the next section "To store the MinION/GridION/PromethION flow cell for later use" before priming and loading the flow cell.</p>	

Flow Cell Wash Kit (EXP-WSH003)

Version: WFC_9088_v1_revE_18Sep2019
Last update: 12/03/2020



Flow Cell Number:

DNA Samples:

INSTRUCTIONS	NOTES/OBSERVATIONS
<input type="checkbox"/> To run a second library straight away, follow the instructions in the "Priming and loading the flow cell" section of the relevant protocol.	
To store the PromethION flow cell for later use	
<input type="checkbox"/> Thaw one tube of Storage Buffer (S) at RT. <input type="checkbox"/> Mix contents thoroughly by pipetting and spin down briefly. <input type="checkbox"/> Rotate the inlet port cover clockwise to reveal the inlet port. A small tract of air may be visible beyond the inlet port. If necessary, using a P1000 draw back a small volume to remove any air (a few µl): <input type="checkbox"/> Set a P1000 pipette to 200 µl <input type="checkbox"/> Insert the tip into the inlet port <input type="checkbox"/> Turn the wheel until the dial shows 220-230 µl, or until you can see a small volume of buffer/liquid entering the pipette tip. <input type="checkbox"/> Slowly add 500 µl of Storage Buffer through the inlet port of the flow cell. <input type="checkbox"/> Close the inlet port cover, and remove any buffer from the waste port at the top of the flow cell. <input type="checkbox"/> The flow cell can now be stored at 4-8° C. <input type="checkbox"/> When you wish to reuse the flow cell, remove from storage, and allow it to warm to RT for ~5 minutes.	